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“MY BEST ASSISTANT ... **CHEMICALLY RESISTANT, PHYSICALLY RESISTANT ...** OR BETTER PUT: UNBEATABLE.”

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I | LABORATORY GLASS BOTTLES AND ACCESSORIES

Due to their high chemical and thermal-shock resistance, DURAN® laboratory bottles are indispensable for daily work in the laboratory. As neutral glass with high hydrolytic resistance, DURAN® belongs to glass type I in accordance with EP, JP and USP.

The extensive range of original accessories includes caps for the widest possible range of applications. DURAN® laboratory bottles are completed by suitable pouring rings from different plastics, which enable drip-free working.

Alongside clear glass laboratory bottles, amber (absorbs light up to approx. 500 nm) and plastic-coated (absorbs light up to approx. 380 nm) bottles are also available to protect light-sensitive substances. The PU plastic coating also provides protection against scratching and, in the event of breakage, retains contents and glass splinters.

Usage tips:

- DURAN® laboratory bottles are, with the exception of the pressure-resistant DURAN® pressure plus bottles, not designed for use under pressure or vacuum conditions.
- When sterilising / autoclaving, the screw cap should

only be loosely applied since with a closed bottle pressure equalisation does not take place. An ideal complement is the membrane cap (see pages 18, 30).

- When cleaning in a dishwasher, load such that the glass bottles, especially the threads, are not in contact with each other.
- The printed scale on many items of DURAN® laboratory glassware is indicative with an accuracy of $\pm 10\%$. Therefore the items are not suitable for use as volumetric glassware.
- DURAN® protect bottles can be autoclaved. The maximum permissible temperature is $+135\text{ °C}$ (max. 30 Min.).
- When freezing media, store the bottles in an inclined position and do not fill more than three quarters full. Bear in mind that plastic caps and pouring rings have restricted service temperature ranges.

DURAN® laboratory glass and reagent bottles are provided with a retrace code. Using the eight-character code and the corresponding article number, a batch and quality certificate for the DURAN® laboratory glass bottles can be obtained at www.duran-group.com.

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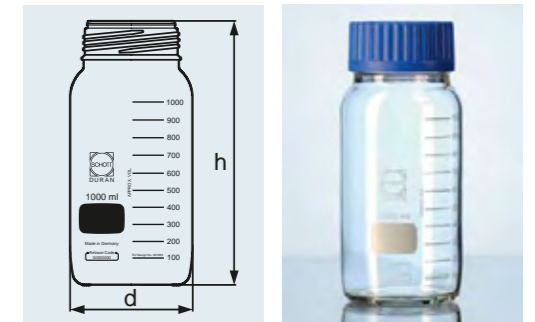
With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. Complete with blue quick release closure (PP¹, integral lip seal) and pouring ring (PP¹) for drip-free pouring and clean, safe working. Service temperature limit of closure and pouring ring: $+140\text{ °C}$. Special thread means opening takes less than a turn. The 80 mm wide outer diameter of the bottle neck permits easy filling and pouring out of powders and viscous substances.

Typical applications: storage, transport, safekeeping and sampling of substances, easy to use with granulated material, powders and viscous media, sampling of hot media.

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DURAN® GLS 80 LABORATORY BOTTLE, WIDE NECK with GLS 80 thread

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DURAN® GLS 80 LABORATORY BOTTLE, WIDE NECK, AMBER
with GLS 80 thread

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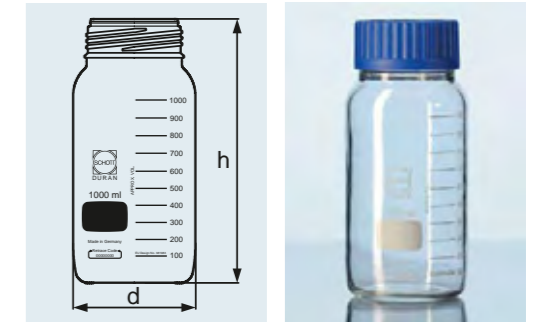
¹ For chemical and thermal resistance, see page 208.

With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. Complete with blue quick release closure (PP¹, integral lip seal) and pouring ring (PP¹) for drip-free pouring and clean, safe working. Service temperature limit of closure and pouring ring: +140 °C. Alongside easy handling, UV protection up to 500 nm. Unchanged DURAN® properties within the bottle, as colouration is only to the outer surface. Very uniform, durable and chemically resistant amber colour due to use of innovative technology.

Typical applications: storage, transport and safekeeping of light-sensitive substances, easy to use with granulated material, powders and viscous media.

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DURAN® GLS 80 PROTECT LABORATORY BOTTLE
with GLS 80 thread, plastic coated

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With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. Complete with blue quick release closure (PP¹, integral lip seal) and pouring ring (PP¹) for drip-free pouring and clean, safe working. Service temperature limit of closure and pouring ring: +140 °C. Service temperature limit of the PU¹ plastic coating: -30 °C to +135 °C. The coating provides scratch, leak and splinter protection and is ideally suited to both the transport and storage of hazardous media or valuable samples. UV protection up to approx. 380 nm wavelength. High transparency. Suitable for microwaving.

Typical applications: storage, transport and safe handling of hazardous substances.

Best.-Nr. Cat. No.	Inhalt Capacity ml	Gewinde Thread GLS	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
11 601 52	500	80	101	148	10
11 601 63	1 000	80	101	218	10
11 601 64	2 000	80	136	248	10
11 601 65	5 000	80	182	310	1

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Permits opening and closing of the DURAN® GLS 80 bottle with only a three-quarter turn. A matching PP¹ pouring ring is also available, permitting clean, drop-free use.

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
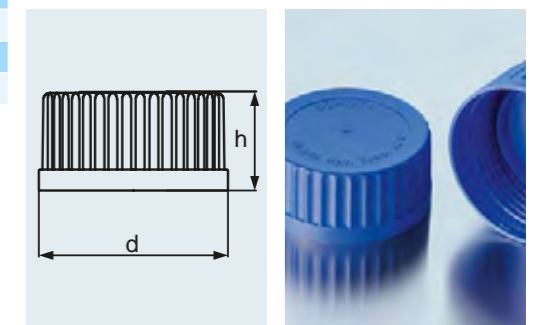
QUICK RELEASE CLOSURE FOR DURAN® GLS 80 LABORATORY BOTTLE
from PP¹, blue, with lip seal

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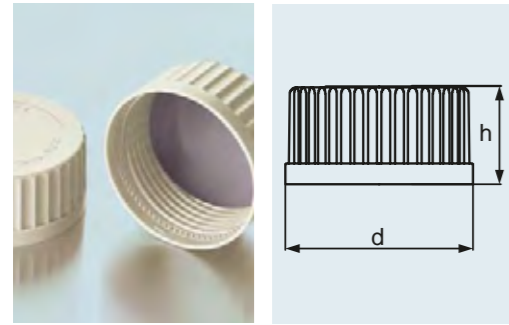QUICK RELEASE CLOSURE FOR DURAN® GLS 80 LABORATORY BOTTLE with sealing disc, (PSU compound¹)

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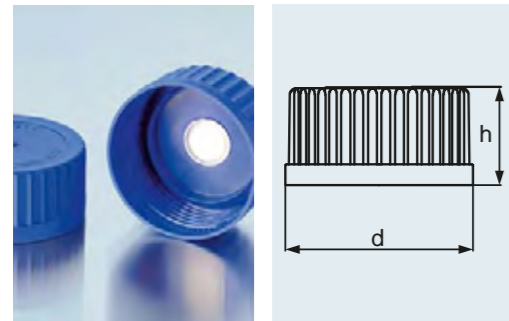
The material used is a special compound based on a polyarylsulphone I. Consequently the chemical, thermal and mechanical properties of the material are noticeably improved and matched to laboratory requirements. Thanks to the thread, the DURAN® GLS 80 bottle can be opened and closed with only a three-quarter turn. The seal, which is coated on both sides with PTFE¹, ensures the bottle can be tightly closed. A matching PTFE¹ pouring ring is also available, permitting clean, drop-free use.

¹ For chemical and thermal resistance, see page 208.

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MEMBRANE SCREW CAP FOR DURAN® GLS 80 LABORATORY BOTTLE from PP¹, blue, with welded-in PTFE¹ membrane for pressure equalisation

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For GLS 80 thread. Ideal for autoclaving processes because the membrane permits pressure equalisation and tight sealing. Hence the risk of contamination is greatly reduced. Ingress of liquids or solids is prevented and the bottle contents remain sterile.

Typical applications: storage or transport of gas generating media, autoclaving of media.

¹ For chemical and thermal resistance, see page 208.

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¹ For chemical and thermal resistance, see page 208.

Materials used: PP¹ and PTFE¹. Flexible modular system. Five different hose diameters (3.2 mm; 6.0 mm; 8.0 mm; 10.0 mm and 12.0 mm) can be connected. Sterile pressure equalisation is possible by using the membrane filter. Unused ports can be provided with a blind cap.

Typical applications: safe transfer of liquid media within a closed and sterile system (evaporation is reduced).

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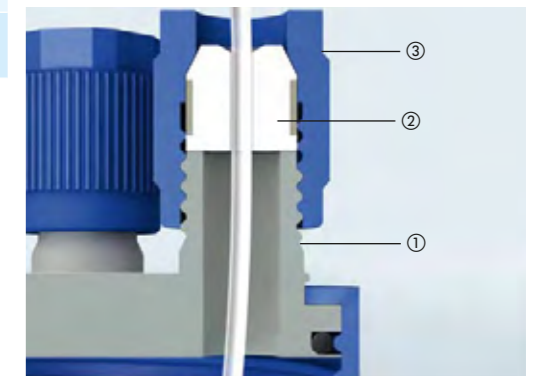
GLS 80 CONNECTION SYSTEM screw cap GLS 80 with four ports (GL 18 thread)

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Schematic diagram of GLS 80 connection system

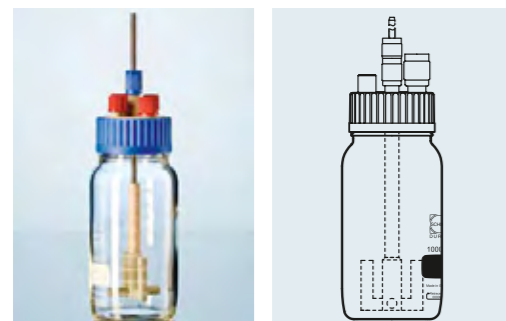
- ① z. B. Best.-Nr. 11 601 76
- ② z. B. Best.-Nr. 11 601 69
- ③ z. B. Best.-Nr. 11 601 74
- ① e.g. Cat. No. 11 601 76
- ② e.g. Cat. No. 11 601 69
- ③ e.g. Cat. No. 11 601 74

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GLS 80 STIRRED REACTOR

materials used: PP¹/PTFE¹/PEEK¹/stainless steel

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The GLS 80 stirred reactor is suitable for a wide range of laboratory mixing processes. The connections (2 x GL 14, 2 x GL 18) provided permit addition or removal of media from the bottle during the mixing process. The whole assembly can be autoclaved and is therefore suitable for use in the biological sector. By using components from the GLS 80 connection system, an additional media bottle (OD hose: 1.6 - 12.0 mm) can be connected or a sterile pressure equalizer attached. Drive for the stirrer is provided by a standard commercial magnetic stirrer. The variable stirrer shaft can be used in DURAN® GLS 80 laboratory glass bottles (1 000 and 2 000 ml) and provides notably improved mixing in comparison with standard magnetic stir bars. The agitating element is exchangeable and use up to 500 rpm is possible.

Typical applications: mixing of liquids, dissolving of solids, simple fermentation processes.

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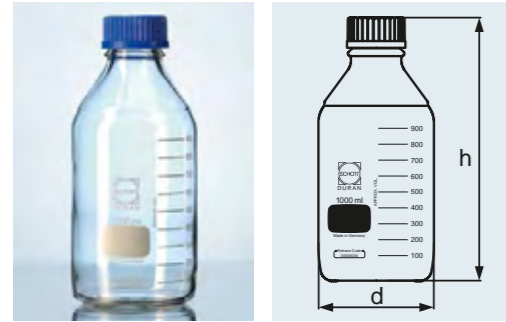
DURAN® LABORATORY BOTTLE
with DIN thread, GL 45

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With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. With proven DURAN® properties. Complete with blue screw cap (PP¹, integral lip seal) and pouring ring (PP¹) for drip-free pouring and clean, safe working. Service temperature level of screw cap and pouring ring: +140 °C. For screw cap assortment, see pages 29–36.

Typical applications: storage, sample preparation, transport.

Best.-Nr. Cat. No.	Inhalt Capacity ml	DIN-Gewinde DIN Thread GL	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
Mit Schraubverschluss und Ausgießring aus PP (blau), klar <i>With screw-cap and pouring ring from PP (blue), clear</i>					
21 801 14 5 ²	25	25	36	74	10
21 801 17 5	50	32	46	91	10
21 801 24 5	100 ³	45	56	105	10
21 801 36 5	250 ³	45	70	143	10
21 801 44 5	500 ³	45	86	181	10
21 801 54 5	1 000 ³	45	101	230	10
21 801 63 5	2 000 ³	45	136	265	10
21 801 73 5	5 000	45	182	335	1
21 801 86 5	10 000	45	227	415	1
21 801 88 5	15 000	45	268	450	1
21 801 91 5	20 000	45	288	510	1
Ohne Verschluss und Ausgießring, klar <i>Without cap and pouring ring, clear</i>					
21 801 14 ²	25	25	36	70	10
21 801 17	50	32	46	87	10
21 801 24	100 ³	45	56	100	10
21 801 36	250 ³	45	70	138	10
21 801 44	500 ³	45	86	176	10
21 801 54	1 000 ³	45	101	225	10
21 801 63	2 000 ³	45	136	260	10
21 801 73	5 000	45	182	330	1
21 801 86	10 000	45	227	410	1
21 801 88	15 000	45	268	445	1
21 801 91	20 000	45	288	505	1

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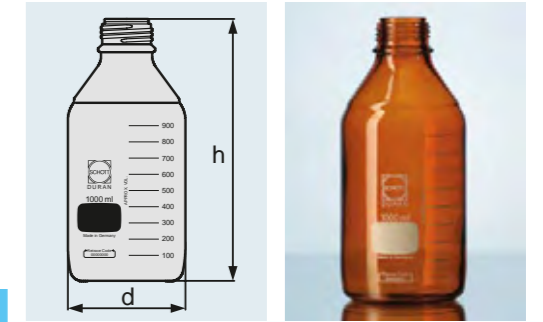
¹ For chemical and thermal resistance, see page 208.

² With specially shaped glass edge for improved pouring out (so that an additional plastic pouring ring is not required).

³ Bottles of 100 – 2 000 ml capacity have a shoulder reinforcement ring corresponding to the bottles nominal capacity.

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DURAN® LABORATORY BOTTLE, AMBER
with DIN thread, GL 45

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With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. UV protection up to approx. 500 nm wavelength. Unchanged DURAN® properties within the bottle, as colouration is only to the outer surface. Very uniform, durable and chemically resistant amber colour due to use of innovative technology. For screw cap assortment, see pages 29–36.

Typical applications: storage and transport of light-sensitive substances.

Best.-Nr. Cat. No.	Inhalt Capacity ml	DIN-Gewinde DIN Thread GL	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
Braun Amber					
21 806 14 ¹	25	25	36	70	10
21 806 17	50	32	46	87	10
21 806 24	100 ²	45	56	100	10
21 806 36	250 ²	45	70	138	10
21 806 44	500 ²	45	86	176	10
21 806 54	1 000 ²	45	101	225	10
21 806 63	2 000 ²	45	136	260	10
21 806 73	5 000	45	182	330	1
21 806 86	10 000	45	227	410	1
21 806 88	15 000	45	268	445	1
21 806 91	20 000	45	288	505	1

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¹ With specially shaped glass edge for improved pouring out (so that an additional plastic pouring ring is not required).

² Bottles of 100 – 2 000 ml capacity have a shoulder reinforcement ring corresponding to the bottles nominal capacity.

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DURAN® PROTECT
LABORATORY BOTTLE
with DIN thread, GL 45, plastic coated

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With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. Service temperature limit of the PU¹ plastic coating: –30 °C to +135 °C (instructions for autoclaving, see page 217). The coating provides scratch, leak and splinter protection and is ideally suited to both the transport and storage of hazardous media or valuable samples. UV protection up to approx. 380 nm wavelength. High transparency. Suitable for microwaving. Caps and pouring rings see pages 29–36.

Typical applications: storage, transport and safe handling of hazardous substances.

Best.-Nr. Cat. No.	Inhalt Capacity ml	DIN-Gewinde DIN Thread GL	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
Ohne Verschluss und Ausgießring, klar Without cap and pouring ring, clear					
10 926 76 ²	25	25	36	70	10
10 926 77	50	32	46	87	10
21 805 24	100 ³	45	56	100	10
21 805 36	250 ³	45	70	138	10
21 805 44	500 ³	45	86	176	10
21 805 54	1 000 ³	45	101	225	10
21 805 63	2 000 ³	45	136	260	10
21 805 73	5 000	45	182	330	1

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DURAN® PRESSURE PLUS
LABORATORY BOTTLE¹
with DIN thread, GL 45

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
With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable ceramic. Pressure resistance conforms to DIN ISO 1595, confirmed by GS marking (TÜV ID: 0000020716). Vacuum and/or pressure resistant from –1 to +1.5 bar due to modified geometry (based on ISO 4796-1). When pressure loaded the following apply: thermal shock resistance 30 K and maximum usage temperature +140 °C. Blue scale for visual differentiation from the standard laboratory bottle. Also available in amber. See pages 29-36 for spare caps and pouring rings.

Typical applications: safe working under pressure or vacuum, sampling under pressure, storage of gas generating media.

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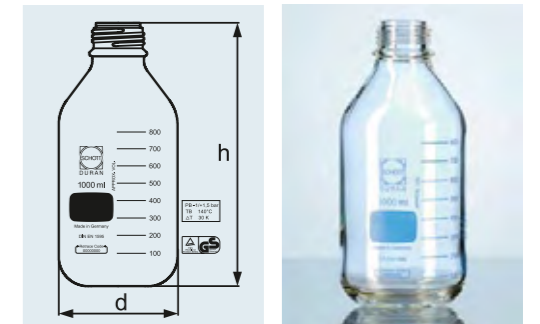DURAN® HPLC BOTTLE
with DIN thread, GL 45

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With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable ceramic. Complete system comprising: DURAN® pressure plus laboratory bottle with 4-port screw cap (PP¹); four screw connections (black, M8 thread) and silicone seals. Connection of different hose diameters (1.6 mm and 3.2 mm) as well as sterile pressure equalisation sets (membrane filter 0.2 µm) is possible. Unused ports can be sealed with silicone blanking seals.

Typical applications: safe transfer of liquid media within a closed and sterile system (evaporation is reduced).

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¹ For chemical and thermal resistance, see page 208.

² Bottle only.

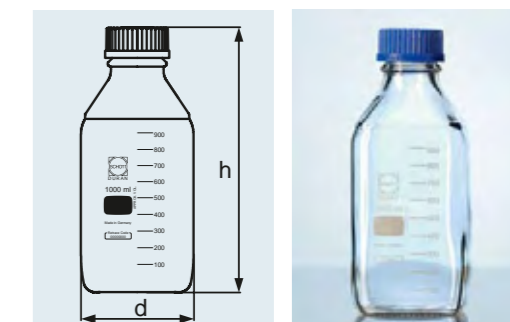
³ Replacement bottle, see page 25.

⁴ Including silicone seals.

⁵ GL 45 screw cap, 4 x M8 screw cap, 12 x silicone seal.

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DURAN® LABORATORY BOTTLE,
SQUARE
with DIN thread, GL 45

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¹ For chemical and thermal resistance, see page 208.

With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. Complete with blue screw cap (PP¹, integral lip seal) and pouring ring (PP¹) for drip-free pouring and clean, safe working. Service temperature limit of screw cap and pouring ring: +140 °C. Ergonomic handling due to angular shape, highly stable, good stackability. Alongside proven DURAN® properties, a space saving of 44% in comparison with standard laboratory bottles (example applies to 100 ml bottles). Screw caps are also available in the following colours: green, yellow and grey (see pages 29–36).

Typical applications: space-saving storage, space-saving transport.

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With easy-to-read scale and large labelling field for easy marking, in fired-on, highly durable white ceramic. Complete with pouring ring and cap from TpCh260¹ (similar to PFA). The premium cap with its PTFE¹ coated silicone seal is colourless and temperature resistant from -196 °C to +260 °C. Together with proven DURAN® properties, TÜV tested thermal shock resistance of 160 K, confirmed by GS-marking (TÜV ID: 0000020715). USP/FDA conformity for the entire system comprised of bottle, screw cap and pouring ring. DMF available upon request (DMF no. 19757). Accurate scale: ± 5%. Additional graduations as well as additional opposing scale simplify reading off.

Typical applications: due to its properties, ideal for applications in the pharmaceutical industry, handling of aggressive media, sterilisation processes (hot air and dry sterilisation) and depyrogenation.

¹ For chemical and thermal resistance, see page 208.

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Compound formulated without colour dye to avoid risk of leaching. Service temperature limit from -196 °C to +260 °C. Very high chemical resistance. High leak tightness through use of PTFE¹ coated silicone seal. DMF (no. 19757, in conjunction with premium bottle) available upon request. A matching TpCh260¹ pouring ring is also available, permitting clean, drop-free use.

Typical applications: due to its properties, ideal for applications in the pharmaceutical industry, handling of aggressive media, demanding sterilisation processes such as hot air or dry sterilisation and depyrogenation.

¹ For chemical and thermal resistance, see page 208.

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The tamper-evident screw cap is available with lip seal or PTFE¹ coated silicone seal. The integral plastic ring tears when the cap is first opened and is retained on the bottle neck. Therefore it can be clearly seen whether the bottle retains its original seal. After first untwisting, can be used as "normal" screw cap.

Typical applications: secure storage and transport/shipping of valuable media.

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
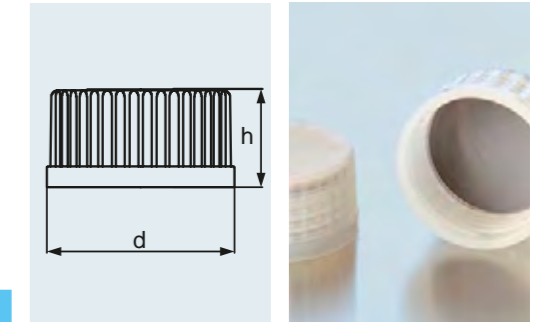
PREMIUM CAP
from TpCh260¹ (similar to PFA) with
PTFE¹ coated silicone seal

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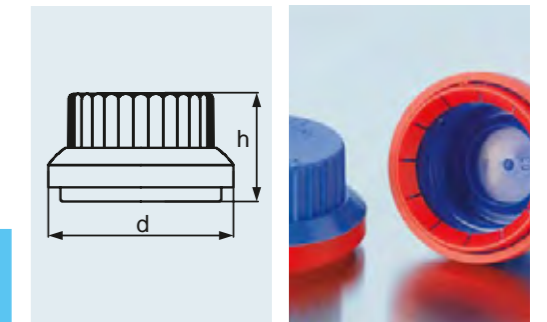
TAMPER-EVIDENT SCREW CAP
from PP¹, for laboratory bottles with DIN thread

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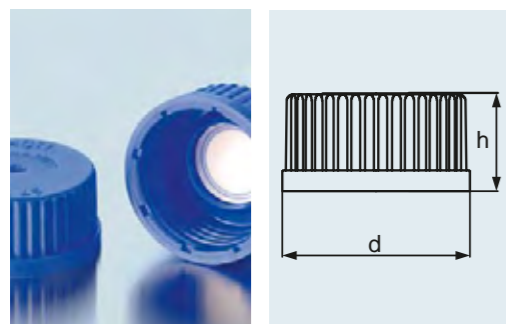MEMBRANE SCREW CAP
from PP¹, blue with welded-in PTFE¹
membrane for pressure equalisation

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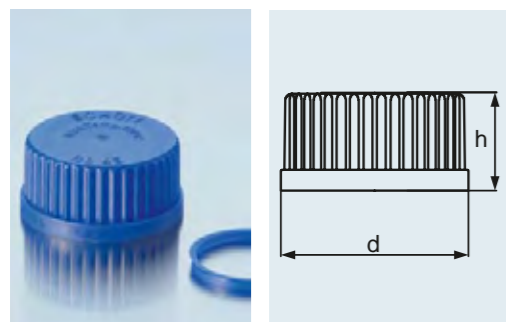
For all GL 45, GL 32 and GL 25 threads.
Ideal for autoclaving processes because
the membrane permits pressure equalisation
and tight sealing, greatly reducing the risk of
contamination. Ingress of liquids or solids is
prevented and the bottle contents remain
sterile.

Typical applications: storage or transport
of gas generating media, autoclaving of
media.

¹ For chemical and thermal resistance, see page 208.

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Available in the colours blue, green, yellow
and grey with matching colour pouring
rings. Distinguishing media types is simplified
and interchanging of screw caps and carry
over of substances is effectively excluded.

Typical applications: visual aid to the safe
identification of different media.

¹ For chemical and thermal resistance, see page 208.

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High leak tightness through use of
PTFE¹ coated silicone seal. More chemically
resistant than PP¹ screw cap. A matching
ETFE¹ pouring ring is also available,
permitting clean, drop-free use.


Best.-Nr. Cat. No.	DIN-Gewinde DIN Thread GL	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
Verschluss Cap ²				
29 240 13	25	33	23	10
29 240 19	32	42	26	10
29 240 28	45	54	28	10
Ausgießring Pouring ring				
29 244 19	32		4	10
29 244 28	45		4	10

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¹ For chemical and thermal resistance, see page 208.

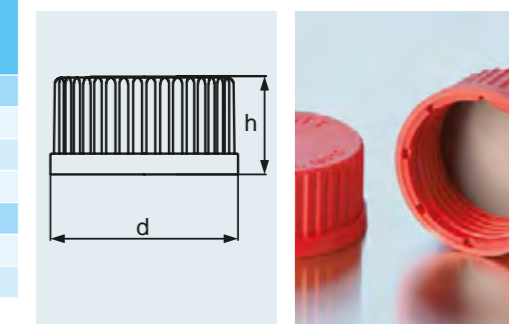² For screw connection caps with aperture, silicone seals
with aperture and septa, see pages 167–169.

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CONNECTION SYSTEM
screw cap GL 45 with two or three ports (GL 14 thread)

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Materials used: PP¹ and PTFE¹. Flexible modular system. Four different hose diameters (1.6 mm; 3.0 mm; 3.2 mm and 6.0 mm) can be connected. Sterile pressure equalisation is possible through use of membrane filter. Unused ports can be provided with a blind cap.

Typical applications: safe transfer of liquid media within a closed and sterile system (evaporation is reduced).

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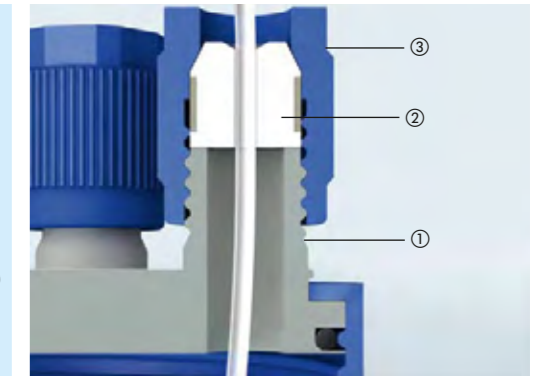
Ordering example

You would like to equip two bottles with a 3-port connection system. You would like to work with two different tube diameters (1.6 mm and 3 mm) and pressure equalization. Consequently, you need the following individual components:

- Screw cap GL 45 3-port (Art. No. 11 297 51) 1 x [2 per pack]
- Screw cap GL 14 (Art. No. 11 298 14) 2 x [2 per pack]
- Insert for screw cap GL14 1.6 mm ID (Art. No. 11 298 15) 2 x [1 per pack]
- Insert for screw cap GL14 3.0 mm ID, (Art. No. 11 298 16) 2 x [1 per pack]
- Pressure equalization set (complete) (Art. No. 11 377 99) 2 x [1 per pack].

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Schematic diagram of GL 45 connection system

- ① z. B. Best.-Nr. 11 297 51
- ② z. B. Best.-Nr. 11 298 15
- ③ z. B. Best.-Nr. 11 298 14

- ① e.g. Cat. No. 11 297 51
- ② e.g. Cat. No. 11 298 15
- ③ e.g. Cat. No. 11 298 14

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DG SAFETY CAPS

screw cap GL 45 with one, two or three ports

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High-quality and flexible connection system, specifically designed for handling aggressive, gas-generating media. Materials used: PP¹, PBT¹ and PTFE¹. All parts coming into contact with media are from PTFE¹. Three different hose diameters (1.6 mm; 2.3 mm; 3.2 mm) can be connected. Pressure equalisation is possible through use of membrane filter. Outgassing of media is reduced by use of a one-way valve that decreases media loss and protects the user from vapours. A free moving PTFE¹ core prevents rotation of the hoses when screwing onto the DURAN[®] bottle. Unused ports can be provided with a blind cap. Optionally available with stopcock to prevent running dry of the hoses or air intakes during bottle changeover. DG safety caps are also available for bottles with standard ground joint.

Typical applications: safe transfer of aggressive media within a closed system (evaporation is reduced).

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The activated carbon filter prevents the escaping of hazardous gases and thus increases safety in the laboratory. Materials used: PBT¹ and PTFE¹. Optional special connection for installations with a leakage system. A free moving PTFE¹ core prevents rotation of the hoses when screwing onto the DURAN[®] bottle. Unused ports can be provided with a blind cap.

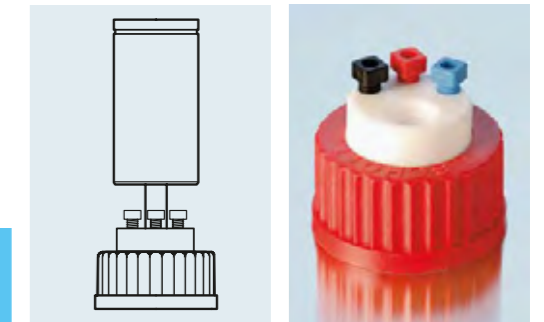
Typical applications: safe storage and collection of liquid laboratory waste.

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DG SAFETY WASTE CAPS

additional connection option for activated carbon filter and leak hose, GL 45

A
121 °C

Tmax.
140 °C

¹ For chemical and thermal resistance, see page 208.

² Complete with GL 45 screw cap, connection screws and conical nipple (ID 3.2 mm, PTFE).

³ Complete with GL 45 screw cap, connection screws, conical nipple (ID 3.2 mm, PTFE) and connection for leak hose.

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¹ For chemical and thermal resistance, see page 208.
² Also suitable for S40 threads.

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Without pouring ring and dust cover cap¹. A maximum usage temperature of +100 °C is recommended. Thermal shock resistance 30 K. Hydrolytic class 3.

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
SCREW CAP BOTTLE, SQUARE, FROM SODA-LIME GLASS narrow neck with thread

Best.-Nr. Cat. No.	Inhalt Capacity ml	Gewinde Thread	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
Klarglas Clear glass					
23 810 24 5	100	32 ²	49	119	10
23 810 36 5	250	32 ²	64	155	10
23 810 44 5	500	32 ²	77	186	10
23 810 54 5	1 000	45 ²	97	223	10
Braunglas Amber glass					
23 816 24 5	100	32 ²	49	119	10
23 816 36 5	250	32 ²	64	155	10
23 816 44 5	500	32 ²	77	186	10
23 816 54 5	1 000	45 ²	97	223	10

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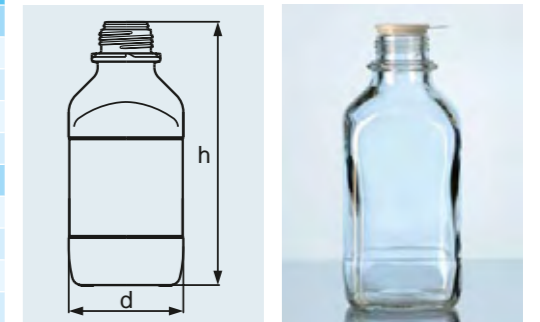
SCREW CAP BOTTLE, SQUARE FROM SODA-LIME GLASS wide neck with thread

Best.-Nr. Cat. No.	Inhalt Capacity ml	Gewinde Thread	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
Klarglas Clear glass					
23 820 17	50	32 ²	48	70	10
23 820 24	100	32 ²	49	111	10
23 820 36	250	45 ²	64	146	10
23 820 44	500	54 ²	76	173	10
23 820 54	1 000	60 ²	97	213	10
Braunglas Amber glass					
23 826 24	100	32 ²	49	111	10
23 826 36	250	45 ²	64	146	10
23 826 44	500	54 ²	76	173	10
23 826 54	1 000	54 ²	97	213	10

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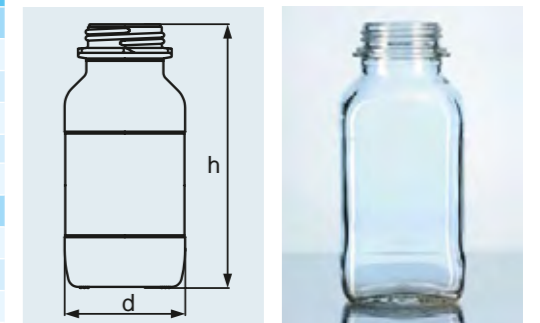SCREW CAP BOTTLE, ROUND, AMBER
FROM SODA-LIME GLASS
with thread

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Complete with pouring ring and dust cover
cap¹. A maximum usage temperature of
+100 °C is recommended. Thermal shock
resistance 30 K. Hydrolytic class 3.


Best.-Nr. Cat. No.	Inhalt Capacity ml	Gewinde Thread	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
Braunglas Amber glass					
23 835 66 5	2 500	45 ²	139	283	1

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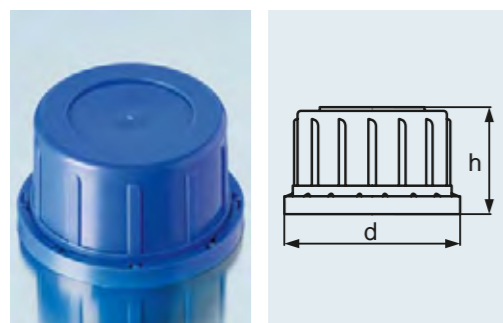TAMPER-EVIDENT SCREW CAP
from PP¹, blue, for soda-lime square screw
cap bottle, narrow neck

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¹ For chemical and thermal resistance, see page 208.
² High form.

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
SECURITY SCREW CAP
from PP¹ with valve², red, for soda-lime
square screw cap bottles, narrow neck

Best.-Nr. Cat. No.	Gewinde Thread	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
29 302 19	32 ³	45	32	10
29 302 28	45 ³	60	35	10

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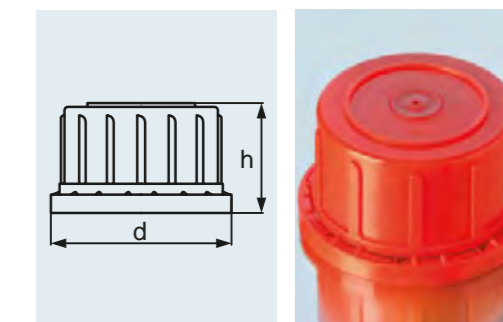¹ For chemical and thermal resistance, see page 208.
² Valve cap opens at 0.1 bar pressure.
³ High form.

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
¹ For chemical and thermal resistance, see page 208.
² Short form.

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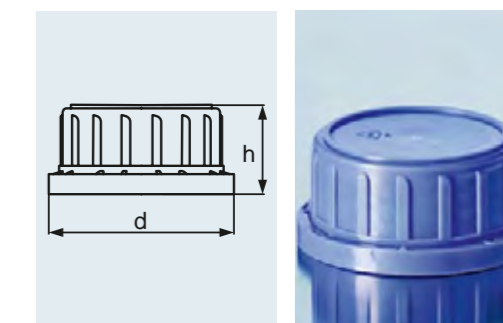TAMPER-EVIDENT SCREW CAP
from PP¹, blue, for soda-lime
square screw cap bottles, wide neck

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DURAN® REAGENT BOTTLE, WIDE NECK
neck with standard ground joint

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DURAN® REAGENT BOTTLE, WIDE NECK
neck with standard ground joint, with standard ground glass flat-head stopper

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¹ No norm available.

With precision ground neck. Only glass components, therefore also suitable for storage of aggressive media, which could attack plastic parts.

Typical application: storage of powders.

With precision ground neck. Only glass components, therefore also suitable for storage of aggressive media, which could attack plastic parts. Also available in amber. Unchanged DURAN® properties within the bottle, as colouration is only to the outer surface. Very uniform amber colour thanks to innovative technology, moreover long durability and high chemical resistance of the amber colouring.

Typical applications: storage of powders.

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With precision ground neck. Only glass components, therefore also suitable for storage of aggressive media, which could attack plastic parts.

Typical application: storage of liquids.

¹ Narrow neck bottles with standard ground joint, 1 000 ml, also suitable for use as burette bottles, see page 113.

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
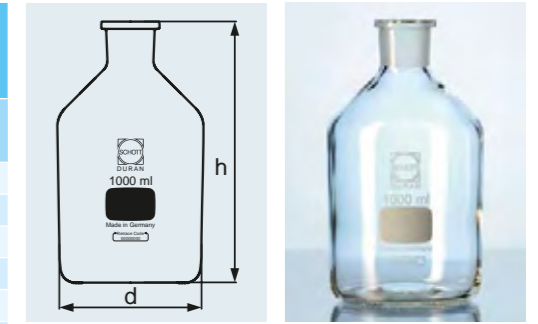
DURAN® REAGENT BOTTLE, NARROW NECK
neck with standard ground joint

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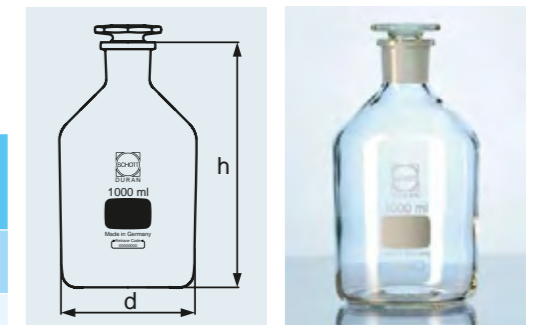
DURAN® REAGENT BOTTLE, NARROW NECK
neck with standard ground joint, with standard ground glass flat-head stopper

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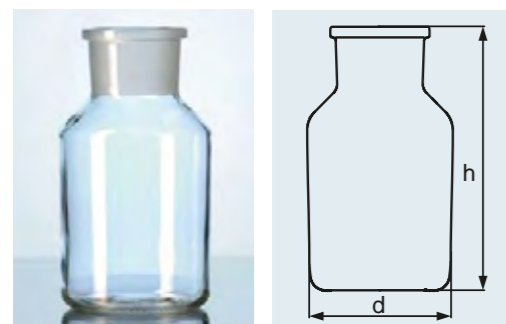REAGENT BOTTLE, WIDE NECK
FROM SODA-LIME GLASS
neck with standard ground joint

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Without stopper¹. A maximum usage temperature of +100 °C is recommended. Thermal shock resistance 30 K. Hydrolytic class 3.

Best.-Nr. Cat. No.	Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	h mm	Hals Neck NS	Verpackungseinheit/ Stück Pack/Quantity
Hals mit Normschliff, ohne Stopfen Neck with standard ground joint, without stoppers						
23 184 17		50	44	79	24/20	10
23 184 24	23 187 24	100	52	97	29/22	10
23 184 36	23 187 36	250	71	129	34/24	10
23 184 44	23 187 44 ²	500	86	164	45/40	10
23 184 54	23 187 54	1 000	107	200	60/46	10

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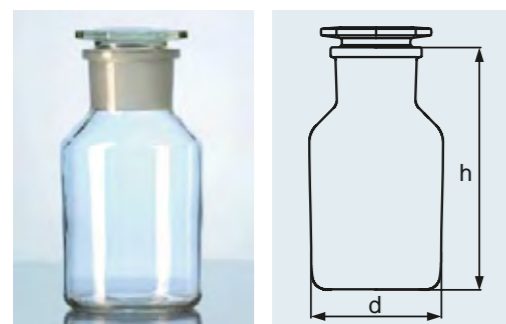
Note on using DURAN® glass stoppers with soda-lime glass bottles: If the bottle and the stopper have a temperature difference greater than 30 °C, the glass stoppers can become stuck!

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REAGENT BOTTLE, WIDE NECK
FROM SODA-LIME GLASS
neck with standard ground joint,
with standard ground glass flat-head stopper¹

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A maximum usage temperature of +100 °C is recommended. Thermal shock resistance 30 K. Hydrolytic class 3.

Best.-Nr. Cat. No.	Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	h mm	Hals Neck NS	Verpackungseinheit/ Stück Pack/Quantity
Hals mit Normschliff und NS-Glasstopfen Neck with standard ground joint and glass flat-head stopper						
23 185 17		50	44	79	24/20	10
23 185 24	23 188 24	100	52	97	29/22	10
23 185 36	23 188 36	250	71	129	34/24	10
23 185 44	23 188 44 ²	500	86	164	45/40	10
23 185 54	23 188 54	1 000	107	200	60/46	10

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¹ Replacement stoppers, see pages 44–45.
² Neck NS 45/27

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REAGENT BOTTLE, NARROW NECK
FROM SODA-LIME GLASS
neck with standard ground joint

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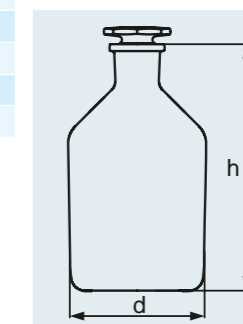
REAGENT BOTTLE, NARROW NECK
FROM SODA-LIME GLASS
neck with standard ground joint,
with standard ground glass flat-head stopper

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GLASS STOPPER, AMBER
with standard ground joint, octagonal

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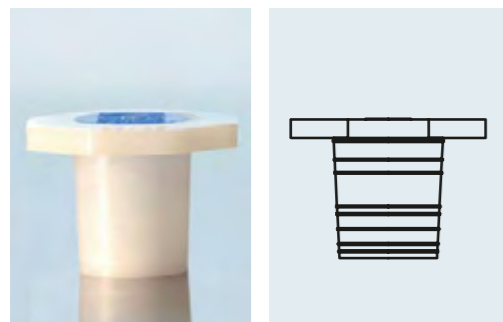
PLASTIC STOPPER
from polyethylene¹

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GLASS STOPPER
with standard ground joint, octagonal

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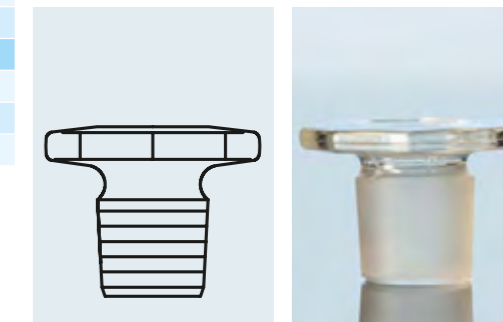
GLASS STOPPER
with short ground joint, octagonal

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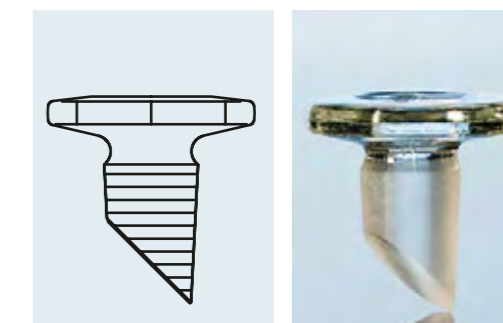
DURAN® GLASS STOPPER
ground conical, for reagent bottles
(oxygen bottles according to Winkler)

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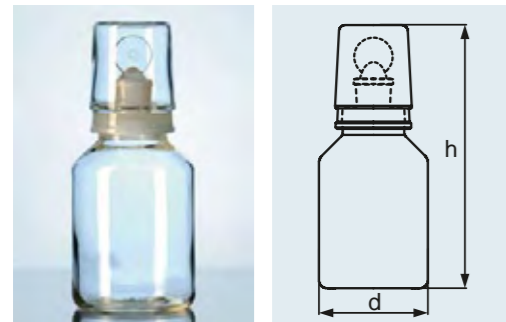
DURAN® LEVELLING BOTTLE
with stopper-bed and olive

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DURAN® ACID BOTTLE¹
with standard ground "pennyhead" stopper,
conical shoulders, interchangeable glass cap

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Dosing of liquids is possible via an outlet.

Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
24 708 36	250	73	130	10
24 708 44	500	89	164	10
24 708 54	1 000	111	200	1

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In addition to the ground stopper, a glass cap with ground joint is supplied. This provides an improved seal and protection against acid vapours.

Best.-Nr. Cat. No.	Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	h mm	NS	Verpackungseinheit/ Stück Pack/Quantity
21 275 24	21 275 24 6	100	55	145	19/17	10
21 275 36	21 275 36 6	250	75	180	19/17	10
21 275 44	21 275 44 6	500	82	220	24/20	10
21 275 54	21 275 54 6	1 000	109	260	29/32	10

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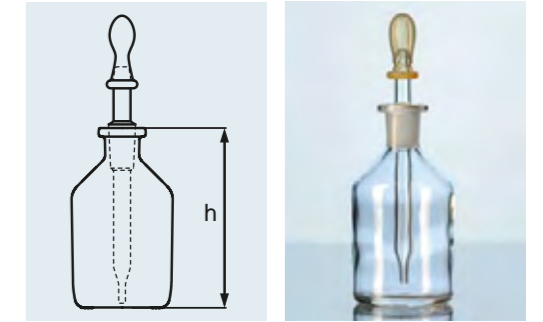
DROPPING BOTTLE
FROM SODA-LIME GLASS
with interchangeable clear glass pipette¹ standard
ground joint, complete with rubber teats²

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
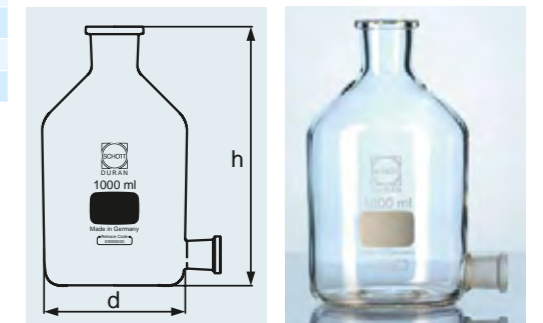
DURAN® ASPIRATOR BOTTLE
tubulated with standard ground joint,
without stoppers, neck unground

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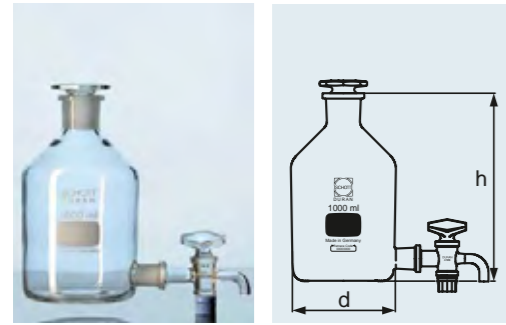DURAN® ASPIRATOR BOTTLE
tubulated with standard ground joint,
complete with standard ground stopcock
and standard ground stopper

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STOPCOCK WITH STANDARD
GROUND JOINT FOR ASPIRATOR
BOTTLE

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DURAN® ASPIRATOR BOTTLE
neck with DIN thread GL 45,
tabulator with GL 32

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Dosing of liquids is possible via a stopcock.

Best.-Nr. Cat. No.	Inhalt Capacity ml	d mm	h mm	Bodentubus Tubulature NS	Hals Neck NS	Orig. Pack./Stück Stand. pack/quantity
24 702 44	500	86	164	19/26	24/29	10
24 702 54	1 000	107	200	19/26	29/32	10
24 702 63	2 000	134	249	19/26	29/32	10
24 702 73	5 000	182	320	29/32	45/40	1
24 702 86	10 000	228	398	29/32	60/46	1
24 702 91	20 000	289	492	29/32	60/46	1

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Complete with screw connection cap¹,
silicone seal and stopcock with PTFE²
spindle. Dosing of liquids is possible via a
stopcock.

Best.-Nr. Cat. No.	Inhalt Capacity ml	DIN-Gewinde DIN Thread GL	d mm	h mm	Verpackungseinheit/ Stück Pack/Quantity
24 703 54	1 000	45	101	225	1
24 703 63	2 000	45	136	260	1
24 703 73	5 000	45	182	330	1
24 703 86	10 000	45	230	410	1

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² For chemical and thermal resistance, see page 208.

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
STOPCOCK FOR ASPIRATOR BOTTLE
with PTFE¹ spindle, for GL 32 screw thread

Best.-Nr. Cat. No.	Bohrung Hole mm	für Inhalt for Capacity ml	Orig. Pack./Stück Stand. pack/quantity
24 147 03	6	1 000 + 2 000	1
24 147 04	8	5 000 + 10 000	1

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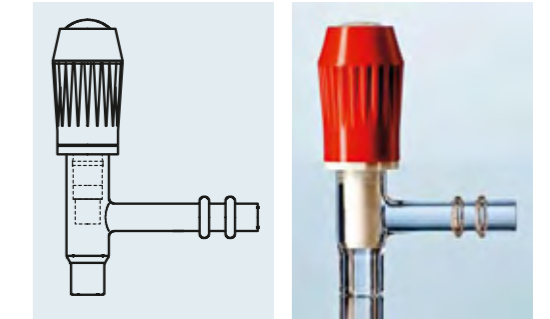PTFE-SILICONE¹ SEAL
for GL 32 screw thread