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“MY LONG-TERM PROJECT ... TO SEPARATE
THE IMPORTANT FROM THE UNIMPORTANT ...
IF ONLY EVERYTHING WAS AS **EASY** AS
SEPARATING USING DURAN®.”

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6 | GLASS FILTRATION APPARATUS AND ACCESSORIES

Due to their high chemical and thermal shock resistance DURAN® filters and the corresponding filter plates are ideal for separations, e. g. with strong acids or alkalis. Thus they offer advantages in comparison with other materials such as plastic or paper. DURAN® filter products have a maximum operating temperature of +450 °C and are therefore far superior to other materials.

The corresponding filtration vessels are specially optimised to the matching filtration apparatus and are vacuum-tight due to their special geometry and high wall thickness. This characteristic has been approved by the German TÜV and marked with the "GS"-indication.

The glass filters are classified as being in porosity classes 0 to 5 according to their nominal maximum pore size. The following table shows the corresponding porosity range. The specified pore sizes always relate to the largest pore in the plate. This specification also characterises the minimum nominal size of particles which may be retained by the filtration.

Usage tips:

- The maximum permissible operating temperature is +450 °C.

- Uniform heating is recommended to avoid thermal stresses and resultant breakages.
- Heat glass filtration apparatus with disk diameters of more than 20 mm in initially cold ovens or sterilisers only.
- The heating or cooling rate should not exceed 8 °C/min.
- When filtering hot substances observe the thermal shock resistance and, if necessary, preheat the filtration apparatus in a drying cabinet.
- Wet filtration apparatus should be heated slowly up to +80 °C and dried for one hour before increasing the temperature further.

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Particular suited to use with hot or aggressive media thanks to the thermal shock and chemical resistance of DURAN®. Also available in **soda-lime glass**.

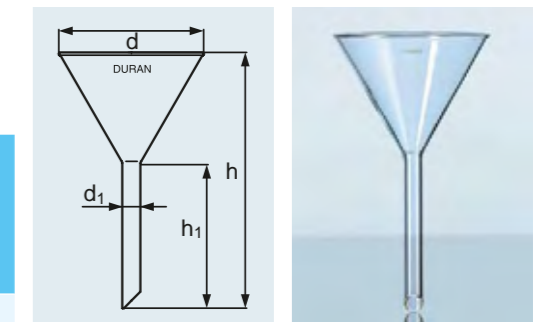
Typical applications: decanting and filtration of substances.

From DURAN® glass with its good thermal shock and chemical resistance. Also available in **soda-lime glass**.

Typical applications: decanting of powdered substances and granulated material.

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
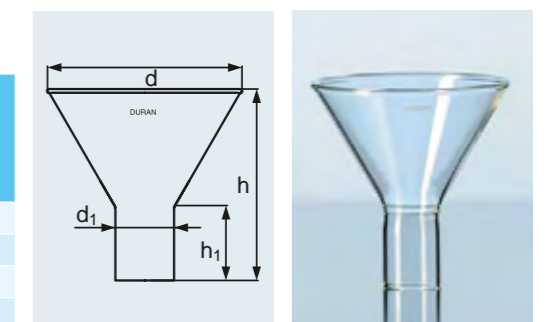
DURAN® POWDER FUNNEL with short, wide stem

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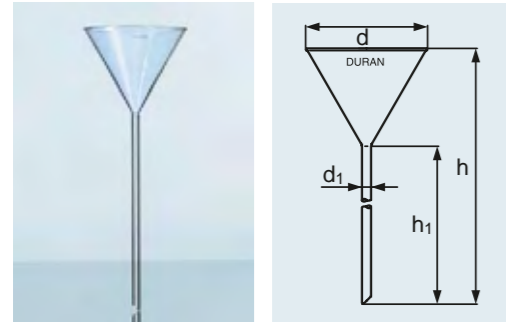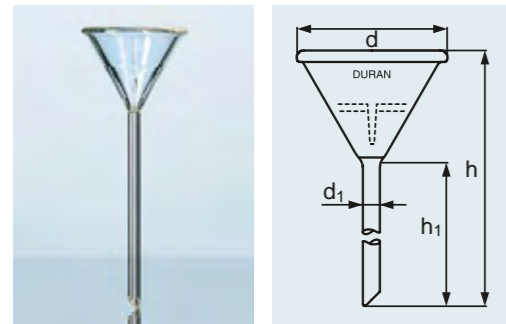
DURAN® ANALYTICAL FUNNEL
for quick filtration

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From DURAN® glass with its good thermal shock and chemical resistance. Also available in soda-lime glass.

Typical applications: filtering and decanting of liquids of different densities.

¹ Non-DIN ISO size.

² Details apply only to DURAN® version.

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From DURAN® glass with its good thermal shock and chemical resistance.

Typical application: for rapid liquid filtration.

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From DURAN® glass with its good thermal shock and chemical resistance. The ribbed form is ideal for filtering with round-paper.

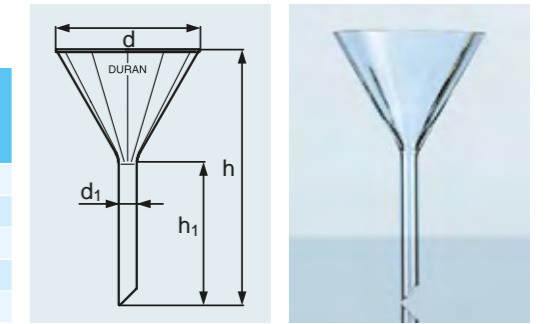
Typical application: filtering of liquids.

Made from DURAN® glass with its good thermal shock and chemical resistance. Filter funnels mate to the filtering flask via a conical rubber seal (GUKO).

Typical application: filtering of liquids.

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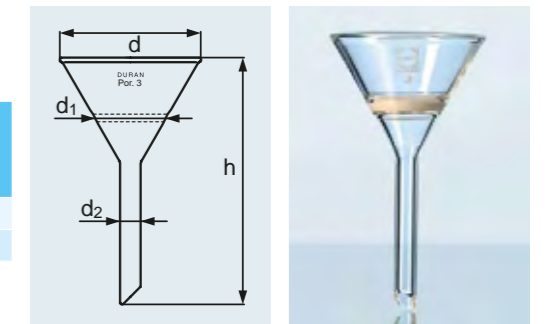
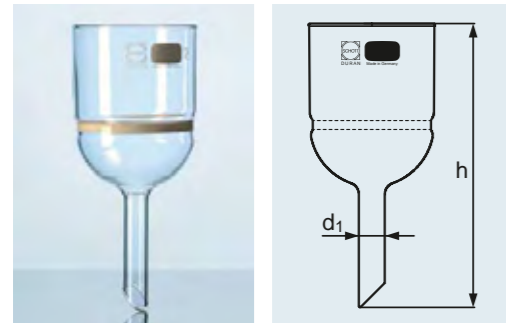
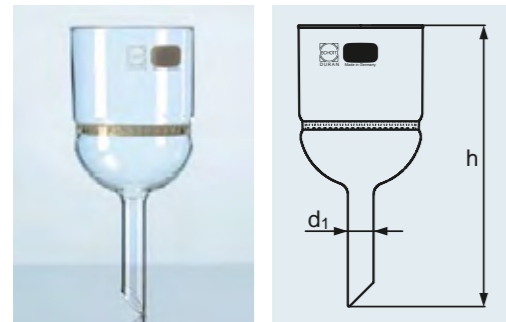
From DURAN® glass with its good thermal shock and chemical resistance. Filter funnels mate to the filtering flask via a conical rubber seal (GUKO).

Typical applications: qualitative inorganic analysis and preparative chemistry.

¹ Substitute required porosity (see page 222) for X in the Cat. No.

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
From DURAN® glass with its good thermal shock and chemical resistance. The Büchner funnel features a glass support for membrane and paper filters.

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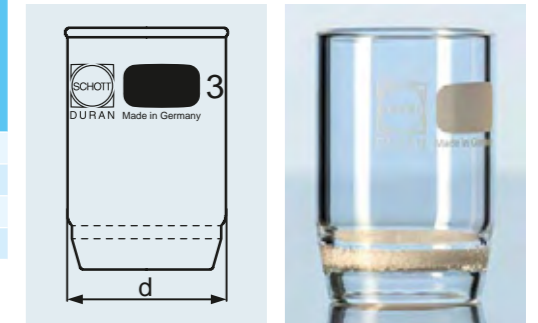
¹ For chemical and thermal resistance, see page 208.

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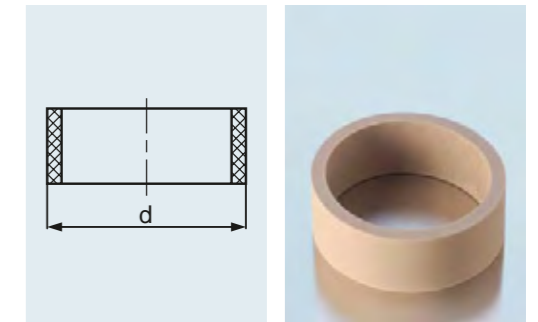
RUBBER SLEEVE
from EPDM¹ for filter crucibles

Best.-Nr. Cat. No.	d mm	d _i mm	h mm	passende Gummi- manschette Best.-Nr. suitable rubber sleeve Cat. No.	Verpackungseinheit/ Stück Pack/Quantity
24 316 16	27	10	108	29 201 14	10
24 316 22	34	10	110	29 201 21	10
24 316 26	41	10	125	29 201 26	10
24 316 32	50	10	132	29 201 31	10


DURAN® FILTRIERVORSTOSS

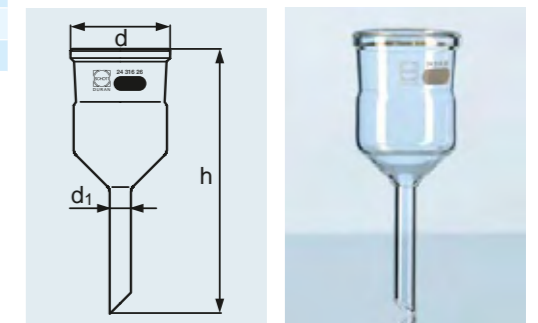DURAN® FILTER CRUCIBLE /
FUNNEL ADAPTER

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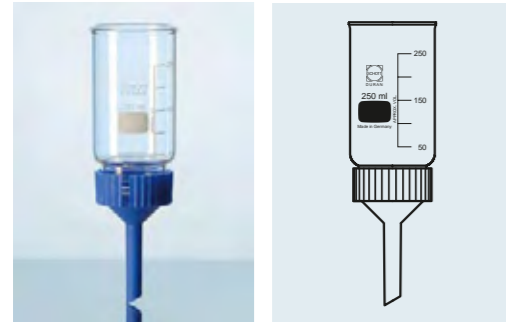DURAN® FILTER FUNNEL HEAD
WITH PP¹ FUNNEL AND TWO
FKM¹ SEALS

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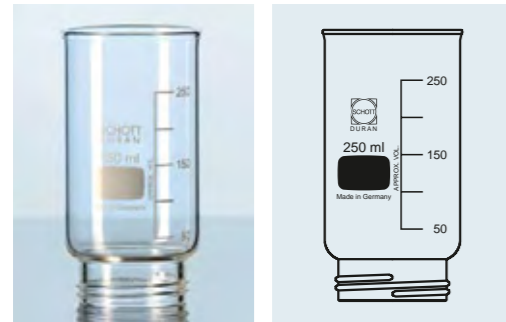
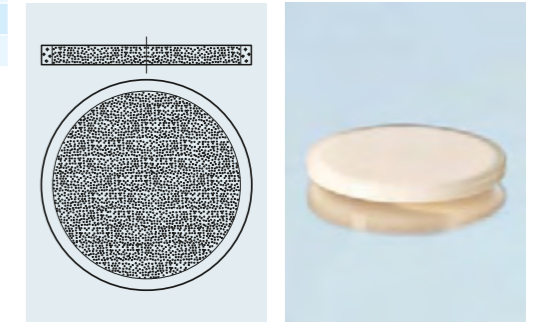
DURAN® FILTER DISK
WITH GLASS RIM
patented

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From DURAN® glass with its good thermal shock and chemical resistance. Fused glass rim.

Best.-Nr. Cat. No.	Porosität Porosity	für Platte for disc d mm	Verpackungseinheit/ Stück Pack/Quantity
25 202 4X ¹	1-4	24	1
25 205 0X ¹	1-4	50	1
25 209 0X ¹	1-4	90	1

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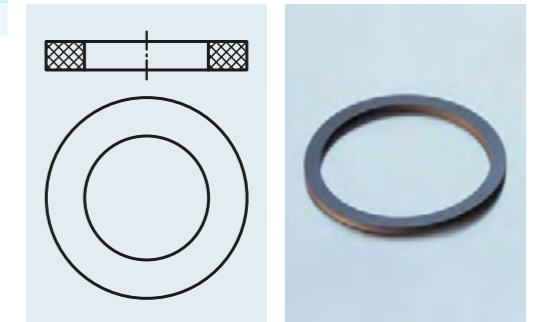
Typical application: in-line filtration of gas lines to remove solid impurities (e.g. dust).


Best.-Nr. Cat. No.	Porosität Porosity	Platte Disc d mm	d ₁ mm	Verpackungseinheit/Stück Pack/Quantity
25 855 0X ¹	1-4	30	14	1
25 855 1X ¹	1-4	60	16	1
25 855 2X ¹	1-4	90	16	1

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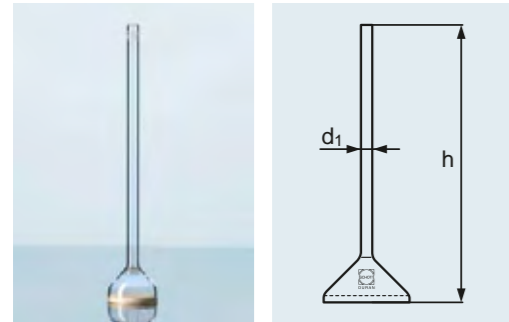DURAN® IMMERSION FILTER
for reverse filtration

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Typical application: extraction of clear filtrate (not the filtrand).


Best.-Nr. Cat. No.	Porosität Porosity	Platte Disc d mm	d ₁ mm	h mm	Verpackungseinheit/Stück Pack/Quantity
25 855 6X ¹	1-4	35	10	210	1

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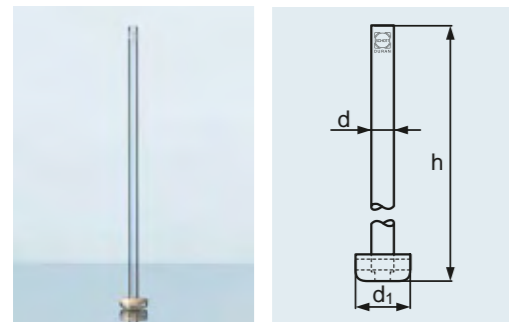Typical application: reactions between gases and liquids.

Best.-Nr. Cat. No.	Porosität Porosity	d mm	d ₁ mm	h mm	Verpackungseinheit/Stück Pack/Quantity
25 856 0X ¹	0-2	6	22	250	5
25 856 1X ¹	0-2	9	25	250	5
25 856 2X ¹	1	11	34	250	5

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With screw-connection system. The insertion height of the head is adjustable. Individual parts can also be ordered separately (see page 128).

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Typical applications: cleaning (“washing”) of gases with solvents.

Best.-Nr. Cat. No.	Inhalt Capacity ml	DIN-Gewinde DIN Thread GL	Olive Hose connection d mm	Verpackungseinheit/ Stück Pack/Quantity
24 713 00	500	45	9	1

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With screw-connection system. The insertion height of the head is adjustable. The distribution of the gas in the liquid is carried out through the filter disk and thus the absorption efficiency significantly increased. Gas washing bottles also work reliably at high flow rates. Individual parts can also be ordered separately (see page 128).

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
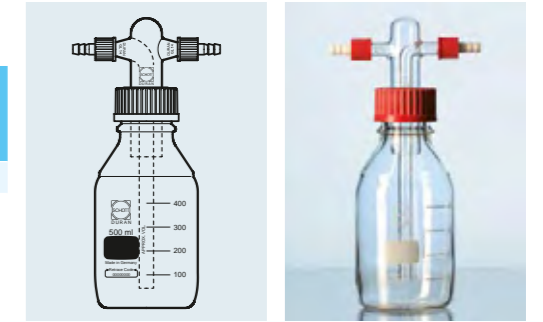
DURAN® GAS WASHING BOTTLE
head Drechsel-type, without filter disk

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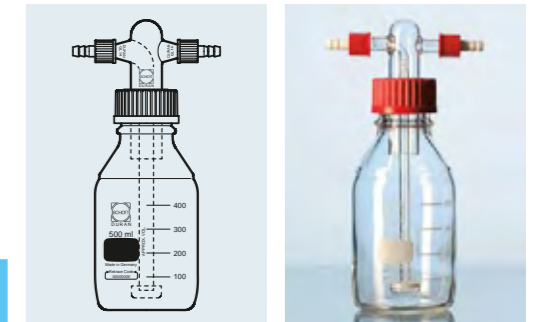DURAN® GAS WASHING BOTTLE
head Drechsel-type, with filter disk

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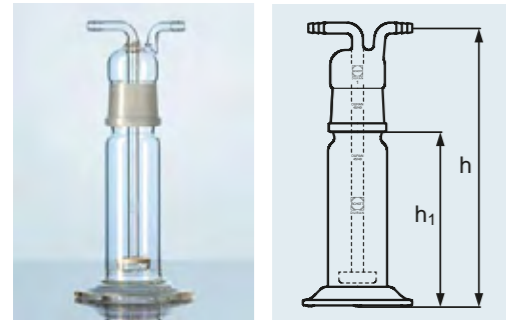
DURAN® GAS WASHING BOTTLE
head with filter disk, with standard ground joint

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Typical applications: cleaning („washing“) of gases with solvents.

Best.-Nr. Cat. No.	Porosität Porosity	Inhalt Capacity ml	Napf Bowl d mm	NS Joint	h mm	h ₁ mm	Olive Hose connection d mm	Verpackungs- einheit/Stück Pack/Quantity
25 701 01	I	100	25	34/35	250	180	10	1
25 702 01	I	250	34	45/40	250	160	10	1

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
Typical application: cleaning („washing“) of gases with solvents.

Best.-Nr. Cat. No.	Porosität Porosity	Inhalt Capacity ml	NS Joint	d mm	h mm	h ₁ mm	Olive Hose connection d mm	Verpackungs- einheit/Stück Pack/Quantity
25 703 01	I	350	29/32	60	250	180	10	1

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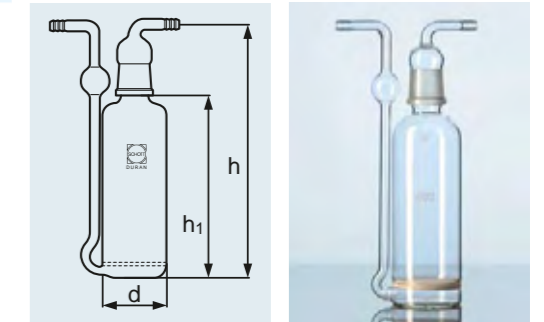DURAN® GAS WASHING BOTTLE
with fused-in filter disk,
with standard ground joint and cap

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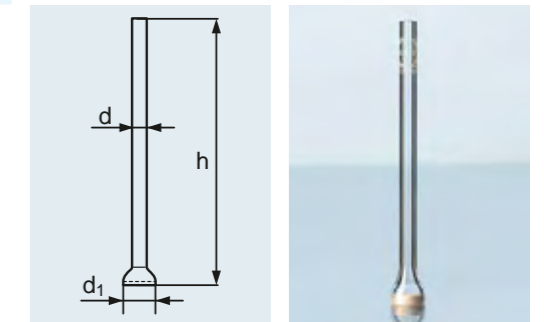
DURAN® MICRO IMMERSION FILTER
for reverse filtration


Best.-Nr. Cat. No.	Porosität Porosity	d mm	h mm	Verpackungseinheit/Stück Pack/Quantity
25 857 2X ¹	I-4	13	25	10

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
DURAN® MIKRO-FILTERKERZE
ohne Rohr

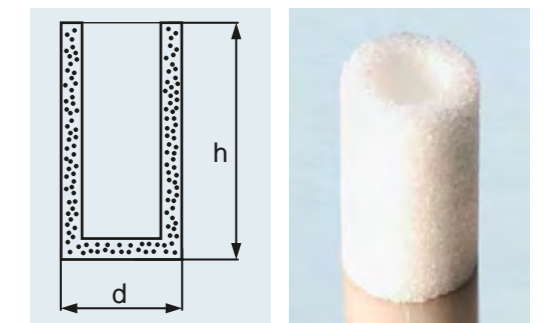DURAN® MICRO FILTER CANDLE
without tube

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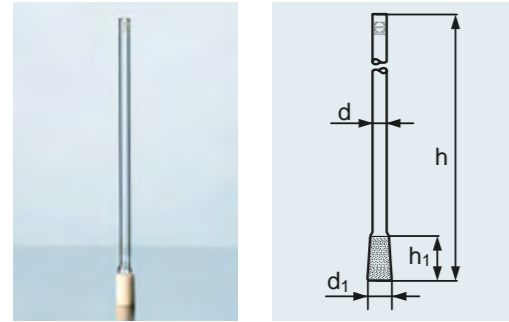DURAN® MICRO FILTER CANDLE
with narrow tube

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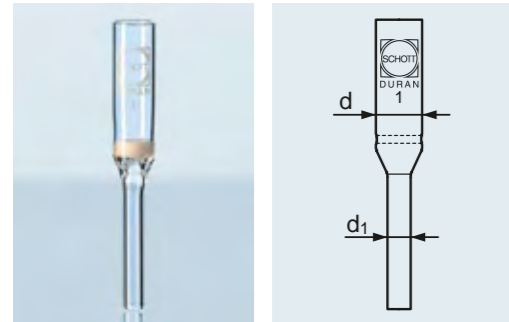
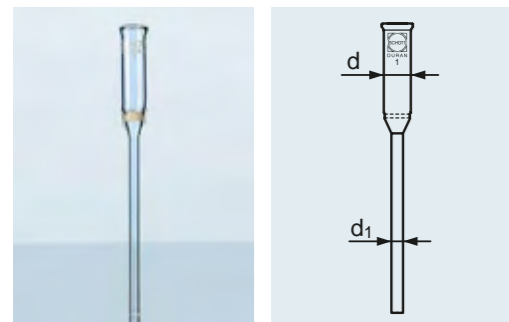
DURAN® MICRO FILTER FUNNEL
Pregl type

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
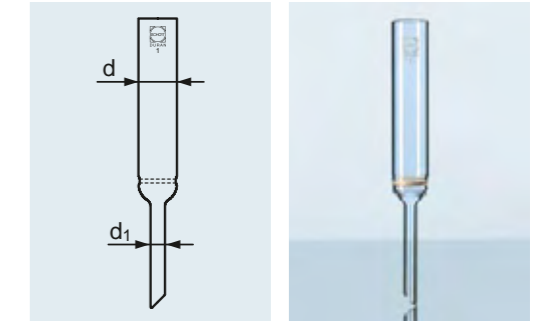
Typical applications: separations by vacuum filtration.

Best.-Nr. Cat. No.	Inhalt Capacity ml	d	d ₁	h	Verpackungseinheit/ Stück Pack/Quantity
21 183 36	250	85	34	155	10
21 183 44	500	105	34	185	10
21 183 54	1 000	135	45	230	10
21 183 63	2 000	166	60	255	1

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Erlenmeyerform

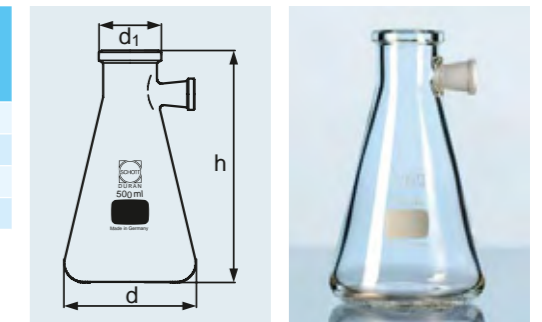
DURAN® FILTERING FLASK
WITH SIDE-ARM SOCKET
Erlenmeyer shape

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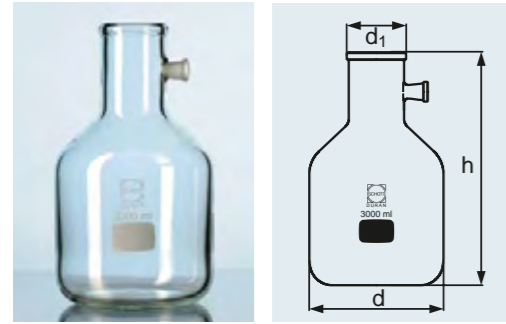DURAN® FILTERING FLASK
WITH SIDE-ARM SOCKET
bottle shape

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Heavy walled for vacuum use. These filtering flasks fulfil the regulations of the "equipment and product safety regulations". For suitable rubber gaskets, see page 135.

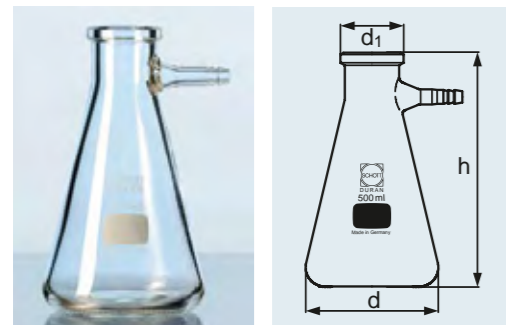
Typical application: separations by vacuum filtration.

Note: These filtering flasks have a ground socket 17.5/26 for vacuum tube of 15 to 18 mm OD (e.g. 6 x 5 mm or 8 x 5 mm, DIN 12 865).

Provision of filtration flasks with a socket has not only made work in preparation and analytical laboratories easier and simpler, but has also reduced the risk of accidents.

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DURAN® FILTERING FLASK
WITH GLASS HOSE CONNECTION
Erlenmeyer shape

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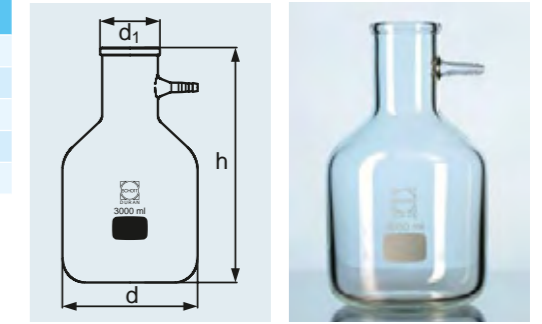
Due to the heavy wall thickness the apparatus is vacuum-tight. For suitable rubber gaskets, see page 135.

Typical application: separations by vacuum filtration.

¹ Does not conform to the "equipment and product safety regulations".

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DURAN® FILTERING FLASK
WITH GLASS HOSE CONNECTION
bottle shape

Heavy walled for vacuum use. For suitable rubber gaskets, see page 135.

Typical application: separations by vacuum filtration.

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
Heavy walled for vacuum use. These filtering flasks fulfil the regulations of the "equipment and product safety regulations". The plastic hose connection is replaceable. For suitable rubber seals, see page 135.

Typical application: Separations by vacuum filtration.

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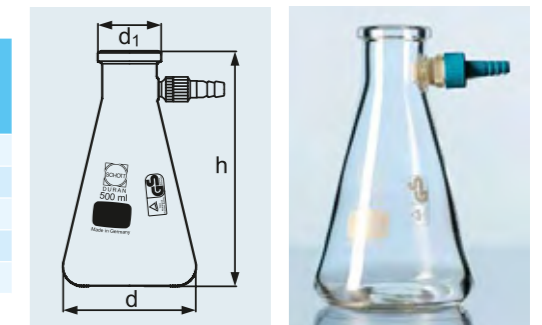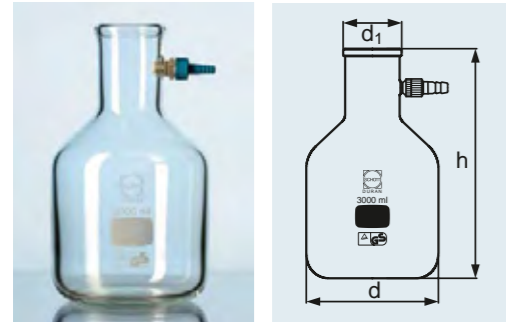
DURAN® FILTERING FLASK
WITH KECK™ ASSEMBLY SET
bottle shape

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
Due to the heavy wall thickness the apparatus is for vacuum use. These filtering flasks fulfil the regulations of the "equipment and product safety regulations". The plastic hose connections can be replaced. For suitable rubber gaskets, see page 135.

Typical application: separations by vacuum filtration.

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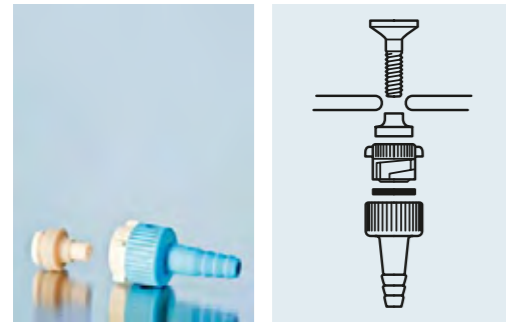With removable plastic hose connection (PBT¹), screw (PP¹), seals (VMQ¹, EPDM¹). Suitable for filtering flasks 100 - 20 000 ml.

Best.-Nr. Cat. No.	Olive Hose connection d mm	Verpackungseinheit/ Stück Pack/Quantity
29 258 54	9	10

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
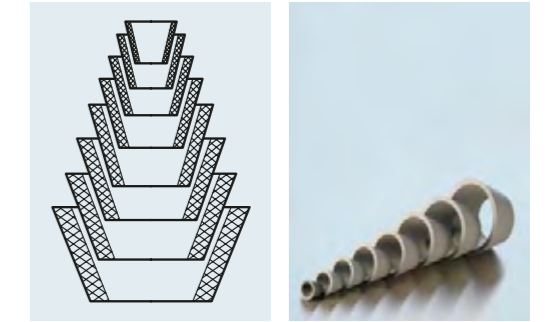
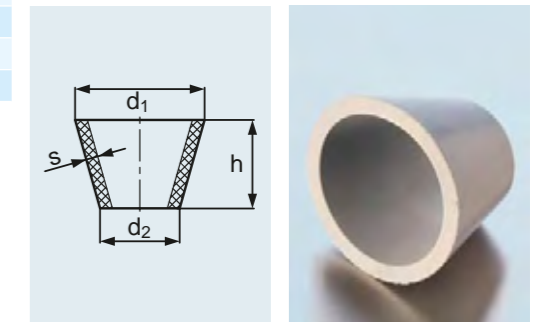
GUKO FROM EPDM¹
conical rubber gaskets for filtering flasks

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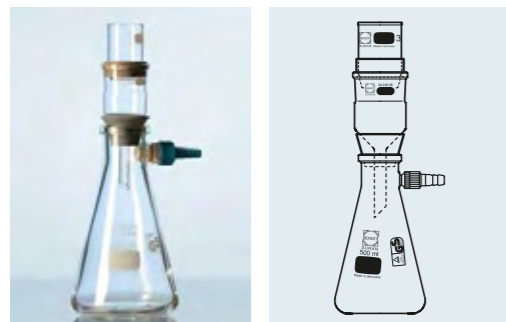DURAN® LID FOR FILTER
APPARATUS, WITT TYPE
with tubulature, standard ground joint 29/32

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¹ Comprises the following order numbers:
 21 204 445, 29 202 27, 24 316 26, 29 201 26,
 25 851 23 and 25 851 24.

Best.-Nr. Cat. No.	DN	Hülse Socket size NS	Verpackungseinheit/ Stück Pack/Quantity
24 398 46	100	29/32	1
24 398 57	150	29/32	1
24 398 61	200	29/32	1

Best.-Nr. Cat. No.	DN	h ₁ mm	h ₂ mm	Hülse Socket size NS	Verpackungseinheit/ Stück Pack/Quantity
24 730 46	100	160	76	29/32	1
24 730 57	150	200	102	29/32	1
24 730 61	200	300	126	29/32	1

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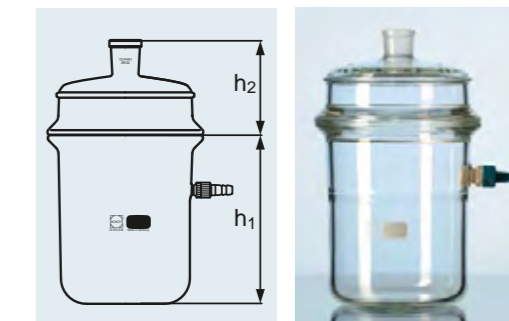
DURAN® FILTER APPARATUS,
WITT TYPE
complete with interchangeable lid and KECK™
assembly set¹, standard ground joint 29/32

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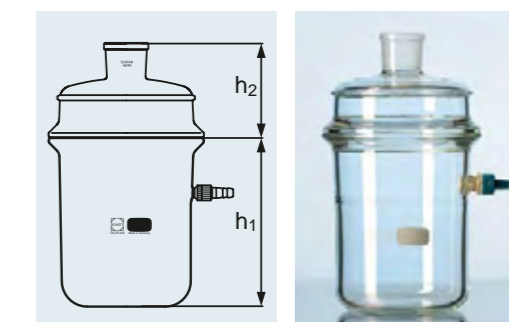
DURAN® FILTER APPARATUS,
WITT TYPE
complete with interchangeable lid and KECK™
assembly set¹, standard ground joint 45/40

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*DURAN® LID FOR FILTER APPARATUS,
 WITT TYPE*
*wide rough-ground tubulature, standard
 ground joint 45/40*

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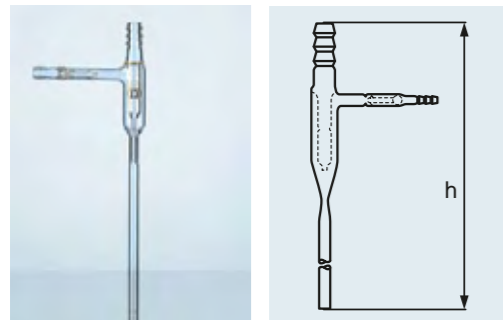
DURAN® WATER JET PUMP
with non-return valve

Best.-Nr. Cat. No.	Wasserverbrauch mind. Water flow rate min. l/h	Wasserdruck mind. Water pressure min. bar	h mm	Verpackungseinheit/ Stück Pack/Quantity
24 362 99	300	1,2	275	1

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*Throughput: 400 l/h at 3.5 bar water
 pressure und 12 °C water temperature.*

Best.-Nr. Cat. No.	Olive passend für Schlauch ID Connection suitable for hose ID mm	Wasserverbrauch mind. Water flow rate min. l/h	Wasserverbrauch mind. Water flow rate min. bar	h mm	Verpackungs- einheit/Stück Pack/Quantity
29 250 01	9-12	170	1	235	1

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WATER JET PUMP
*from plastic (PP¹), with non-return valve,
 hose connection and adapters for 1/2" and 3/4"*